

Kinetic / VLE set-up Intermediate pressure

Purpose

Acid gases can be removed from industrial gas streams with a chemical and/or physical solvent. In this experimental set-up both the reaction kinetics as the acid gas solubility (VLE) can be determined for different gas treating processes. The absorption process will take place in a stainless steel Büchi reactor, which is connected with a high intensity stirrer (gas and liquid phase). The reactor is double walled and connected with a heating bath.


Acid gases

- carbon dioxide, hydrogen sulphide, carbonyl sulphide, mercaptans (methyl- upto pentylmercaptans)

Solvents

- (alkanol)amines, alkaline salts, amino acid salts, ammonia, physical solvents

Picture

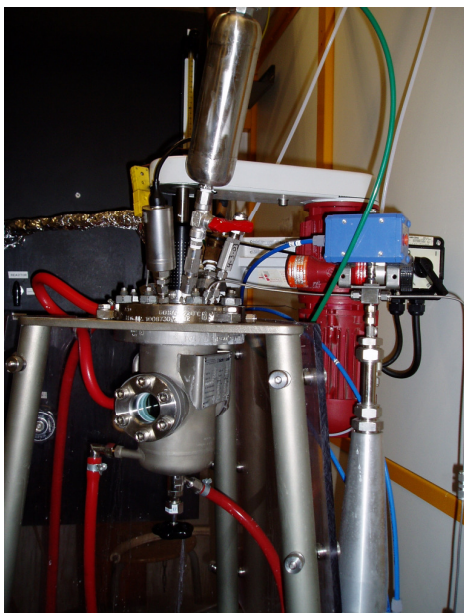

Specifications

100 mbar < P < 35 bar
5 < T < 120 °C
Charge quantity: ± 500 ml

Results

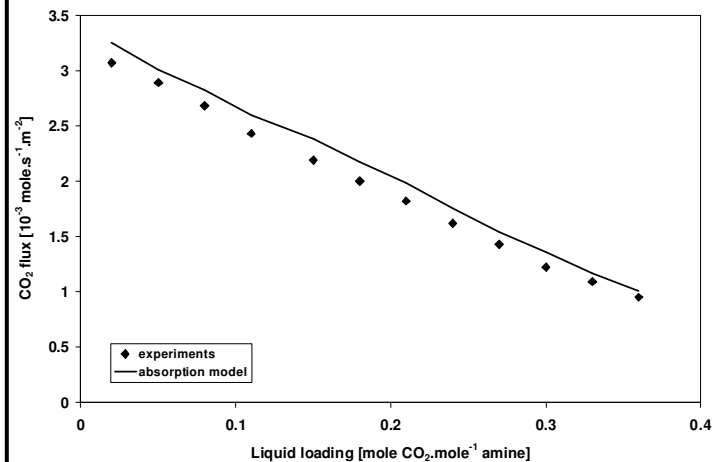


Figure: CO₂ absorption rate in 2 M aqueous DEA at 293 K at different liquid loadings (ref. 1)

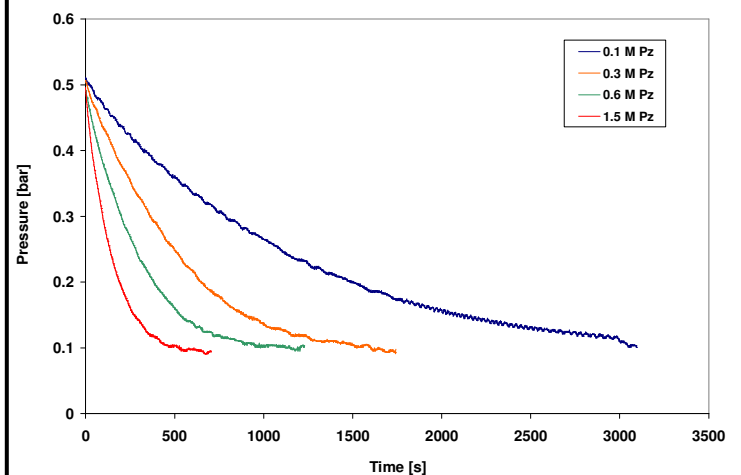


Figure: COS absorption in aqueous piperazine at 313 K (ref.2)

Ref.1 A. Mohan, E. van Elk, S. van Loo, G.F. Versteeg, CHISA 2008, Prague, Czech Republic

Ref.2 P.J.G. Huttenhuis, A. Mohan, S. van Loo, G.F. Versteeg Distillation and Absorption Conference 2006, London, UK

Contact

P. Derks + 31 53 7112 513

P. Huttenhuis + 31 53 7112 508